## CONTACT SLIDE

ITALIANO

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## CONTACT SLIDE

ENGLISH

Flex slide for surfaces bacteriological control  
(ISO 18593:2004)

| CONTACT SLIDE 1                                                  |                                                            |                                                                |        |       |
|------------------------------------------------------------------|------------------------------------------------------------|----------------------------------------------------------------|--------|-------|
| DESCRIPTION                                                      | MEDIA                                                      |                                                                | REF.   | TESTS |
| Slide for total bacterial count and Enterobacteriaceae detection | PLATE COUNT AGAR + TTC + NEUTRALIZING<br>(Bacterial count) | V.R.B.G. AGAR + NEUTRALIZING<br>(Enterobacteriaceae detection) | 525262 | 20    |
|                                                                  |                                                            |                                                                | 53526  | 120   |

| CONTACT SLIDE 2                                                 |                                                            |                                                                     |        |       |
|-----------------------------------------------------------------|------------------------------------------------------------|---------------------------------------------------------------------|--------|-------|
| DESCRIPTION                                                     | MEDIA                                                      |                                                                     | REF.   | TESTS |
| Slide for total bacterial count and yeasts and moulds detection | PLATE COUNT AGAR + TTC + NEUTRALIZING<br>(Bacterial count) | ROSE BENGAL CAF AGAR + NEUTRALIZING<br>(Yeast and moulds detection) | 525272 | 20    |
|                                                                 |                                                            |                                                                     | 53527  | 120   |

| CONTACT SLIDE 3                                             |                                                            |                                                 |        |       |
|-------------------------------------------------------------|------------------------------------------------------------|-------------------------------------------------|--------|-------|
| DESCRIPTION                                                 | MEDIA                                                      |                                                 | REF.   | TESTS |
| Slide for total bacterial count and Staphylococci detection | PLATE COUNT AGAR + TTC + NEUTRALIZING<br>(Bacterial count) | VOGEL JOHNSON AGAR<br>(Staphylococci detection) | 525302 | 20    |
|                                                             |                                                            |                                                 | 53530  | 120   |

| CONTACT SLIDE 4                                            |                                                                 |                                                                      |        |       |
|------------------------------------------------------------|-----------------------------------------------------------------|----------------------------------------------------------------------|--------|-------|
| DESCRIPTION                                                | MEDIA                                                           |                                                                      | REF.   | TESTS |
| Slide for <i>Pseudomonas</i> , yeasts and moulds detection | CETRIMIDE AGAR+ NEUTRALIZING<br>( <i>Pseudomonas</i> detection) | ROSE BENGAL CAF AGAR + NEUTRALIZING<br>(Yeasts and moulds detection) | 525282 | 20    |
|                                                            |                                                                 |                                                                      | 53528  | 120   |

| CONTACT SLIDE 5                                                |                                                                |                                                                     |        |       |
|----------------------------------------------------------------|----------------------------------------------------------------|---------------------------------------------------------------------|--------|-------|
| DESCRIPTION                                                    | MEDIA                                                          |                                                                     | REF.   | TESTS |
| Slide for Enterobacteriaceae and faecal Streptococci detection | V.R.B.G. AGAR + NEUTRALIZING<br>(Enterobacteriaceae detection) | BILE ESCULIN AGAR + NEUTRALIZING<br>(Faecal Streptococci detection) | 525312 | 20    |
|                                                                |                                                                |                                                                     | 53531  | 120   |

| CONTACT SLIDE 8                                        |                                                            |                                                      |        |       |
|--------------------------------------------------------|------------------------------------------------------------|------------------------------------------------------|--------|-------|
| DESCRIPTION                                            | MEDIA                                                      |                                                      | REF.   | TESTS |
| Slide for total bacterial count and coliform detection | PLATE COUNT AGAR + TTC + NEUTRALIZING<br>(Bacterial count) | V.R.B.L. AGAR + NEUTRALIZING<br>(Coliform detection) | 525622 | 20    |
|                                                        |                                                            |                                                      | 53562  | 120   |

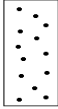
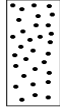
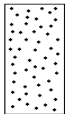
| CONTACT SLIDE 9                                                    |                                                            |                                                            |        |       |
|--------------------------------------------------------------------|------------------------------------------------------------|------------------------------------------------------------|--------|-------|
| DESCRIPTION                                                        | MEDIA                                                      |                                                            | REF.   | TESTS |
| Slide for total bacterial count with inactivation of disinfectants | PLATE COUNT AGAR + TTC + NEUTRALIZING<br>(Bacterial count) | PLATE COUNT AGAR + TTC + NEUTRALIZING<br>(Bacterial count) | 525322 | 20    |
|                                                                    |                                                            |                                                            | 53532  | 120   |

### INSTRUCTIONS FOR USE

1. Unscrew the cap containing the slide. Avoid any contact with the agar surfaces.
2. Flex the cap forming a 90° angle and push for 10 seconds the contact slide on the surface to examine.
3. Rescrew the slide into its tube and incubate at  $36 \pm 1^\circ\text{C}$  for 24-48 hours. For yeasts detection, extend the incubation of further 24-48 hours at  $25-30^\circ\text{C}$ .
4. After incubation examine the bacterial growth on media and estimate the bacterial concentration using the provided scheme.

The interpretation of hygiene level of the surface depends on the examined environment. The following example is referred to work surfaces directly in touch with foods.

#### HYGIENE LEVEL

| Level                  | VERY GOOD                                                                         | GOOD                                                                              | SATISFACTORY                                                                      | DOUBTFUL                                                                            | UNSATISFACTORY                                                                      |
|------------------------|-----------------------------------------------------------------------------------|-----------------------------------------------------------------------------------|-----------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|
|                        |  |  |  |  |  |
| CFU/side               | 1                                                                                 | 2-4                                                                               | 5-13                                                                              | 14-42                                                                               | >42                                                                                 |
| CFU/100cm <sup>2</sup> | 8                                                                                 | 16-32                                                                             | 40-104                                                                            | 112-336                                                                             | >336                                                                                |
| * CFU/10m <sup>2</sup> | 12500                                                                             | 25000-50000                                                                       | 62500-162500                                                                      | 175000-525000                                                                       | >525000                                                                             |

\* 1 CFU/side = 1 x 1.25 x 10<sup>4</sup> CFU/10m<sup>2</sup>

#### CULTURAL RESPONSE

| MEDIUM                        |                     | <i>Escherichia coli</i> | <i>Proteus spp</i> | <i>Klebsiella spp</i> | <i>Pseudomonas spp</i> | <i>Streptococcus spp</i> | <i>Staphylococcus spp</i> | Yeasts/Moulds |
|-------------------------------|---------------------|-------------------------|--------------------|-----------------------|------------------------|--------------------------|---------------------------|---------------|
| <b>V.R.B.G. AGAR</b>          | Colonies appearance | RED-PURPLE              | RED-PURPLE         | RED-PURPLE            | COLORLESS              | INHIBITION               | INHIBITION                | INHIBITION    |
|                               | Medium colour       | RED-PURPLE              | RED-PURPLE         | RED-PURPLE            | RED                    |                          |                           |               |
| <b>V.R.B.L. AGAR</b>          | Colonies appearance | RED-PURPLE              | COLORLESS          | RED-PURPLE            | COLORLESS              | INHIBITION               | INHIBITION                |               |
|                               | Medium colour       | RED-PURPLE              | COLORLESS          | RED-PURPLE            | RED                    |                          |                           |               |
| <b>ROSE BENGAL CAF AGAR</b>   | Colonies appearance | INHIBITION              | INHIBITION         | INHIBITION            | PARTIAL INHIBITION     | INHIBITION               | INHIBITION                | PINK -WHITE   |
|                               | Medium colour       |                         |                    |                       |                        |                          |                           | PINK          |
| <b>VOGEL JOHNSON AGAR</b>     | Colonies appearance | INHIBITION              | PARTIAL INHIBITION | INHIBITION            | INHIBITION             | INHIBITION               | BLACK                     |               |
|                               | Medium colour       |                         |                    |                       |                        |                          | YELLOW                    |               |
| <b>BILE ESCULIN AGAR</b>      | Colonies appearance | INHIBITION              | PARTIAL INHIBITION | PARTIAL INHIBITION    | INHIBITION             | WHITE                    | INHIBITION                |               |
|                               | Medium colour       |                         |                    |                       |                        | BLACK                    |                           |               |
| <b>CETRIMIDE AGAR</b>         | Colonies appearance |                         |                    |                       | GREEN                  |                          |                           |               |
|                               | Medium colour       |                         |                    |                       | BIANCO                 |                          |                           |               |
| <b>PLATE COUNT AGAR + TTC</b> | Colonies appearance | RED                     | RED                | RED                   | RED                    | RED                      | RED                       | RED           |
|                               | Medium colour       | YELLOW                  | YELLOW             | YELLOW                | YELLOW                 | YELLOW                   | YELLOW                    | YELLOW        |

#### STORAGE

10-25°C

#### SHELF LIFE

9 months

#### REFERENCES

- ISO 18593:2004. Microbiology of food and animal feeding stuffs- horizontal methods for sampling techniques from surfaces using contact plate and swabs. ISO/TC 34/SC 9 N 374, march, 1999.
- Handbook of Microbiological Media. Second Edition.1997
- APHA (American Public Health Association) 1970.

#### TABLE OF SYMBOLS

|            |                  |                                                                                     |                         |                                                                                     |                                   |                                                                                       |                                         |
|------------|------------------|-------------------------------------------------------------------------------------|-------------------------|-------------------------------------------------------------------------------------|-----------------------------------|---------------------------------------------------------------------------------------|-----------------------------------------|
| <b>LOT</b> | Batch code       |  | Do not reuse            |  | Manufacturer                      |  | Use by                                  |
| <b>REF</b> | Catalogue number |  | Temperature limitations |  | Contains sufficient for <n> tests |  | Caution, consult accompanying documents |

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**LIOFILCHEM® S.r.l.**